



and 4 ml of 0.02%  $\text{Na}_2\text{S}_2\text{O}_3$   
Incubate for 10 minutes

8. Wash ddH<sub>2</sub>O 2 times 2 minutes

9. Stop 50%MeOH, 12% AcOH Incubate for 10 minutes

10. Wash 50 % MeOH For ~20 minutes at 4°C

Gels may be stored a while at 4°C in 50 % MeOH; or else they may be prepared for drying as follows:

11. Wash 30 % MeOH  
For 30 minutes at 4°C

12. Wash 3% glycerol  
For 30 minutes at 4°C

13. Dry for 1 hour at 80°C between sheets of porous plastic film.